

High Resolution Separations Using Longer Columns and Elevated Temperature

Brian A. Jones, Jody Clark, W. Dale Felix, Stephanie J. Marin

Selerity Technologies, Inc.
2484 W. Custer Road
Salt Lake City, UT 84104

www.selerity.com



High Resolution in HPLC

- Why is it desired
- How it can be accomplished
 - Long columns
 - Small particles
- Ability to move the mobile phase through the column is limiting
 - Raise pressure
 - Decrease viscosity



Column Pressure Drop

The column pressure drop ΔP depends on the mobile phase viscosity η , the mobile phase velocity u (flow rate), column length L and the square of the particle diameter of the packing material d_p :

$$\Delta P = 1000 \eta \cdot u \cdot L / d_p^2$$

$$u = \Delta P \cdot d_p^2 / 1000 \eta \cdot L$$

Increasing plate number in HPLC is thus primarily limited by pressure drop (both L increase and d_p decrease lead to increased pressure drop)




Temperature as a Powerful Tool in LC


- to increase productivity - speed
- for new unique selectivities
- for higher efficiencies (lower mobile phase viscosity)
- for improved peak symmetry
- for higher sensitivities (less band-broadening)
- to reduce organic solvent usage – “green chromatography”



Viscosity of Water vs. Temperature

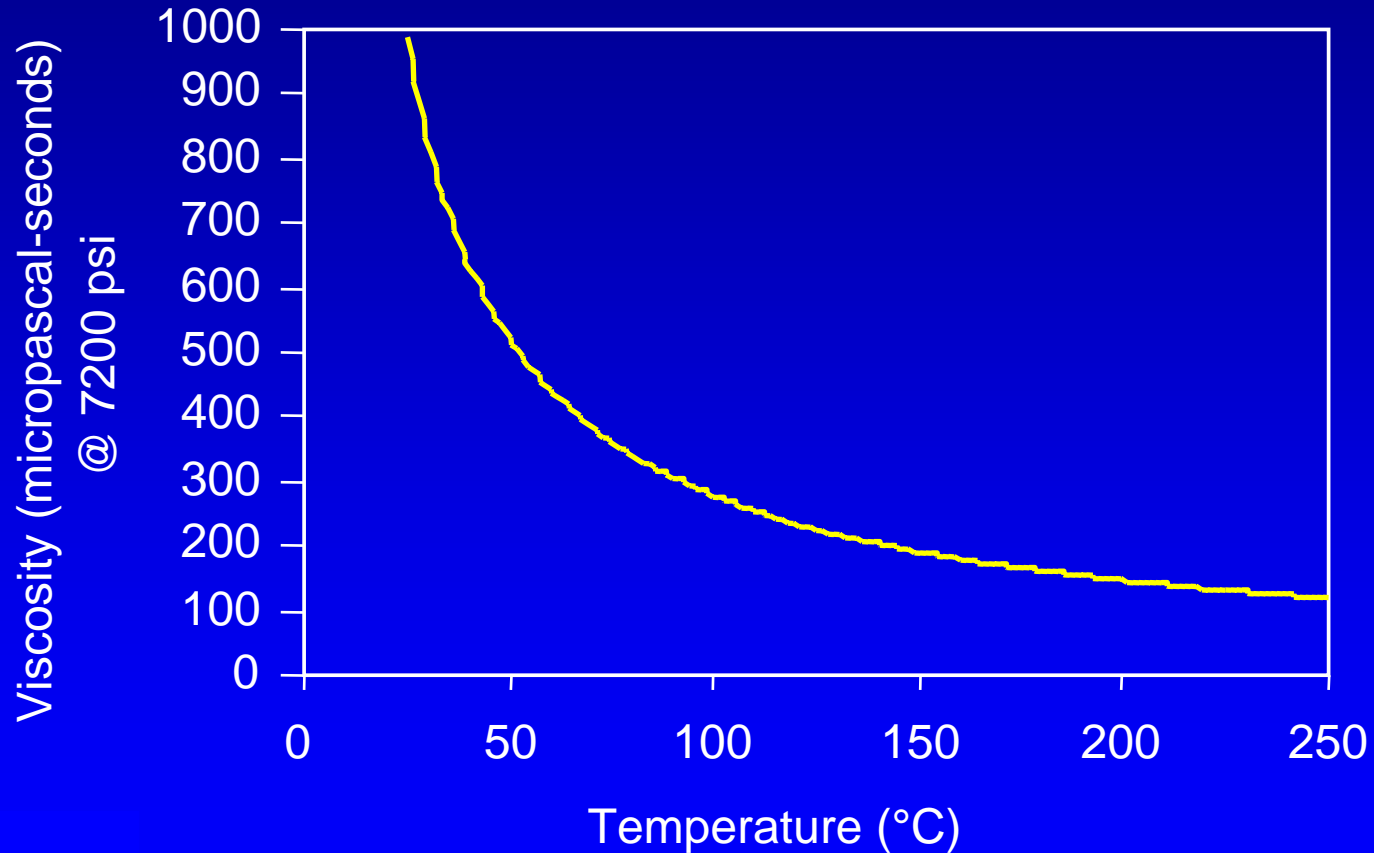

Viscosity Reduction at Elevated Temperature Enables the Use of Long Columns

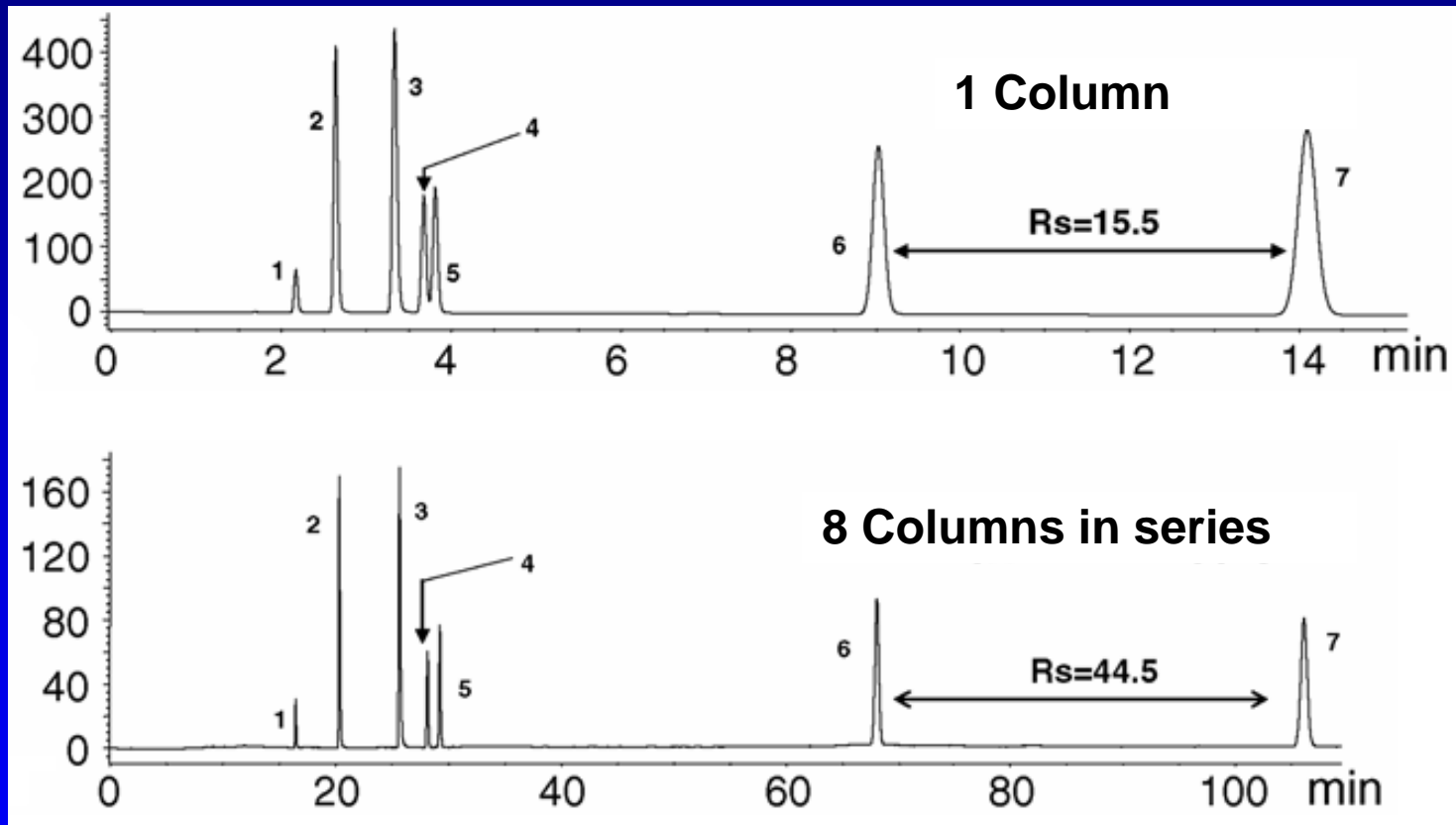

25 cm columns, 5 μ m particles; 80°C; flow rate 1.0 mL/min, 40%ACN. Peaks: 1. uracil; 2. caffeine; 3. pyridine; 4. phenol; 5. aniline; 6. benzene; 7. toluene



Narrow-Bore Columns

- Sensitivity
- Small samples
- Low mobile phase usage
- Detector compatibility
- Fast thermal program tracking




Historical Microbore Columns

- Generally used for normal phase separations only when at elevated temp or moderate pH with reversed-phase solvents
- Packed fused silica capillaries
- PEEK encapsulated or PEEK sealing components (limits temperature to $\sim 100^{\circ}\text{C}$)




Silica Solubility in Water vs. pH

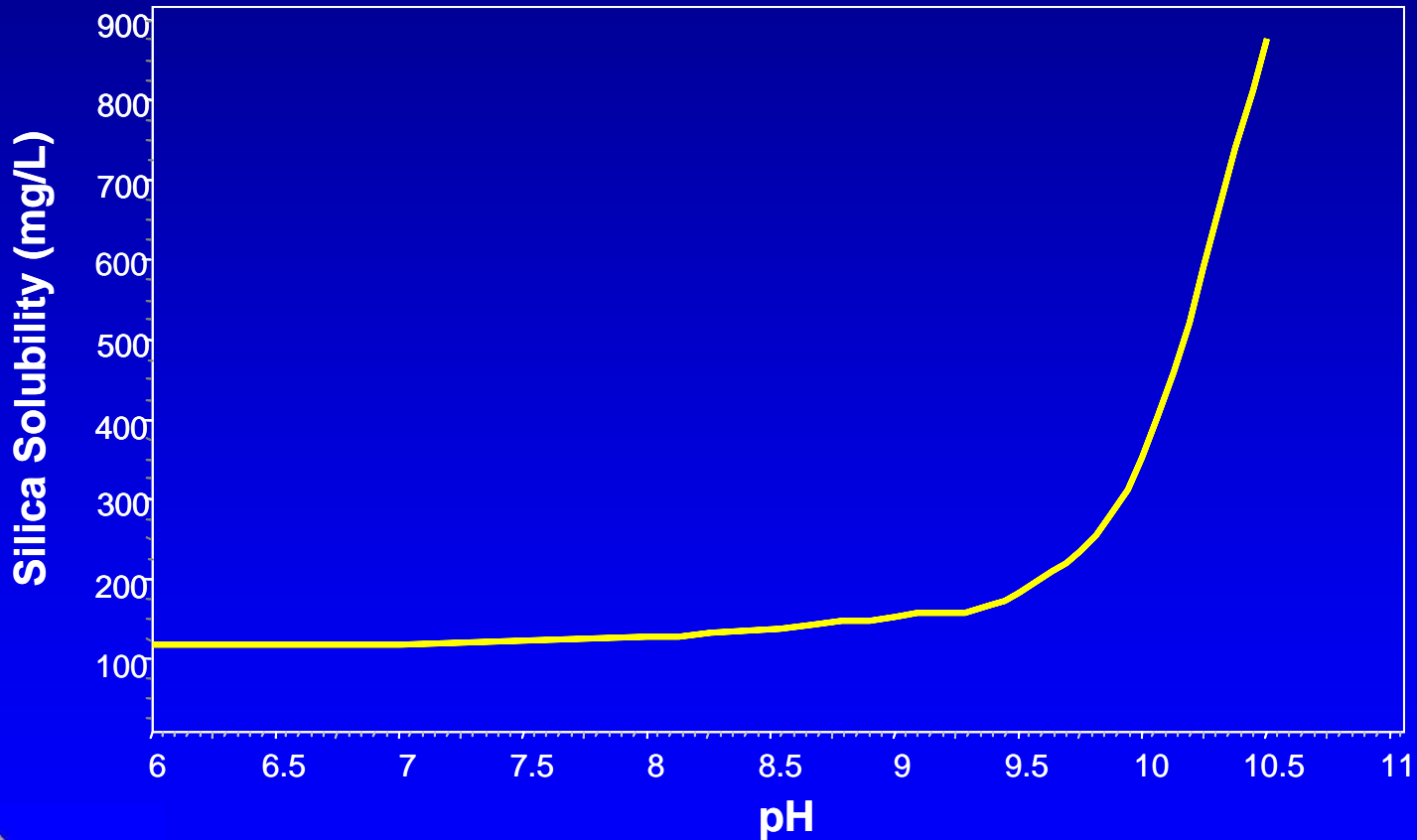

--From R.K. Iler, The Chemistry of Silica, John Wiley & Sons, New York, NY, (1979)

Silica Solubility in Water vs. Temperature

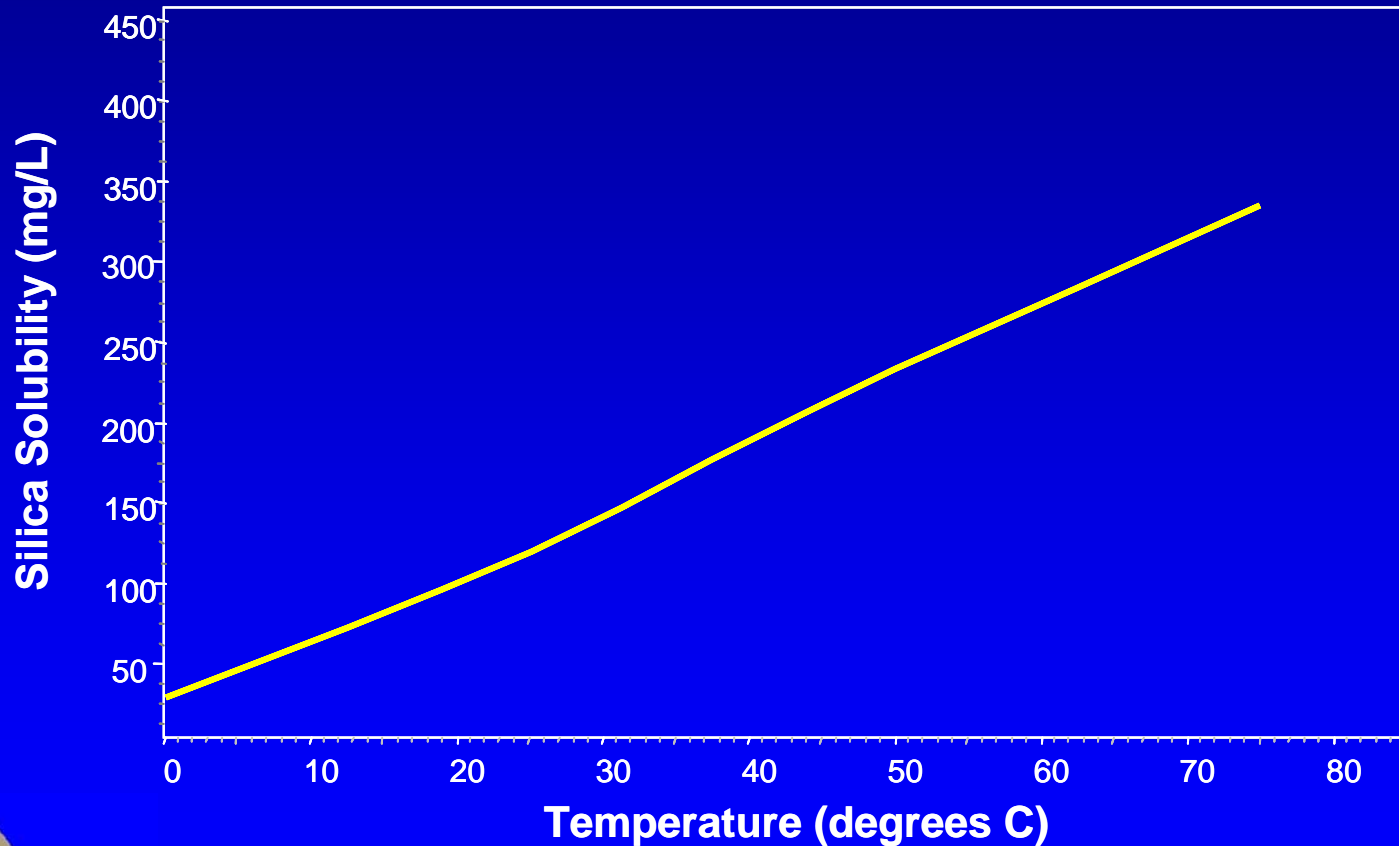

--From R.K. Iler, The Chemistry of Silica, John Wiley & Sons, New York, NY, (1979)

Silica Capillary Dissolution Rate

- More than just bulk solubility in the fluid
- Affected by fluid flow across the surface
- Water concentration and pH influence corrosivity



Comments on the Use of Fused Silica Tubing Based Packed Capillaries for Reversed-Phase Separations Under Aggressive Conditions:

“A silica saturator is used to minimize the dissolution of silica from the capillary wall (not the column, the packing is not silica-based) by the superhot water.”

“Overheating the pre-saturator can result in dissolved silica subsequently depositing in the injector.”

-T. S. Kephart and P. K. Dasgupta, *Talanta* **56** (2002) 977-987



...Bursting of fused silica capillaries occurred even with a silica saturator.

For long term studies, stainless steel capillaries were used...


•T. S. Kephart and P. K. Dasgupta, *Talanta* **56** (2002) 977-987



Selerity SFC Columns



ASTM-5186 Column Performance

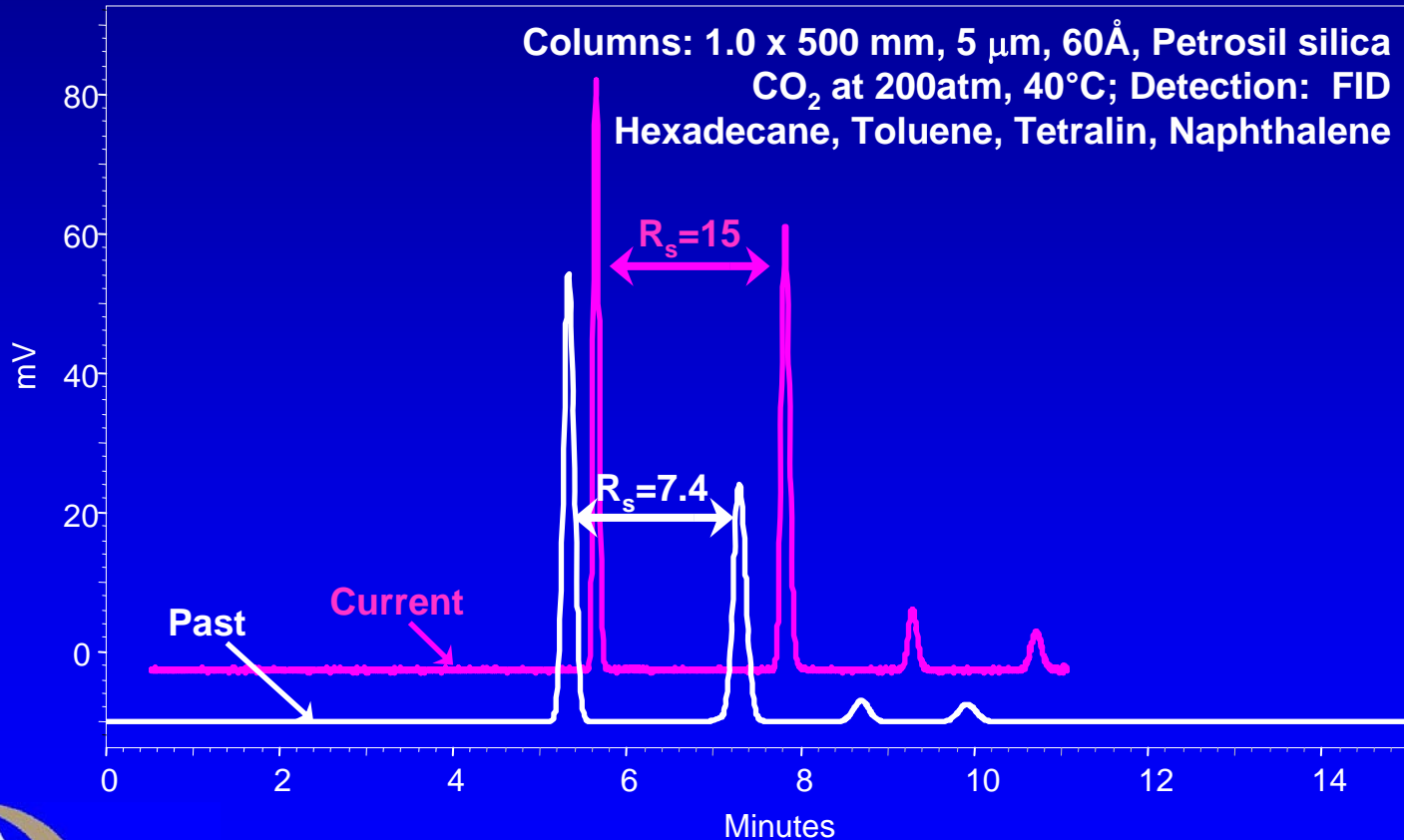


Keys to Improving Narrow-bore Column Performance

- Improvements in packing technology
- Improvements in producing long lengths of highly polished narrow bore stainless steel tubing
- Development of strong particles with traditional silica selectivity and stability in aqueous mobile phases over wide pH and temperature ranges




pHidelity™C₁₈/Blaze₂₀₀™ Surface

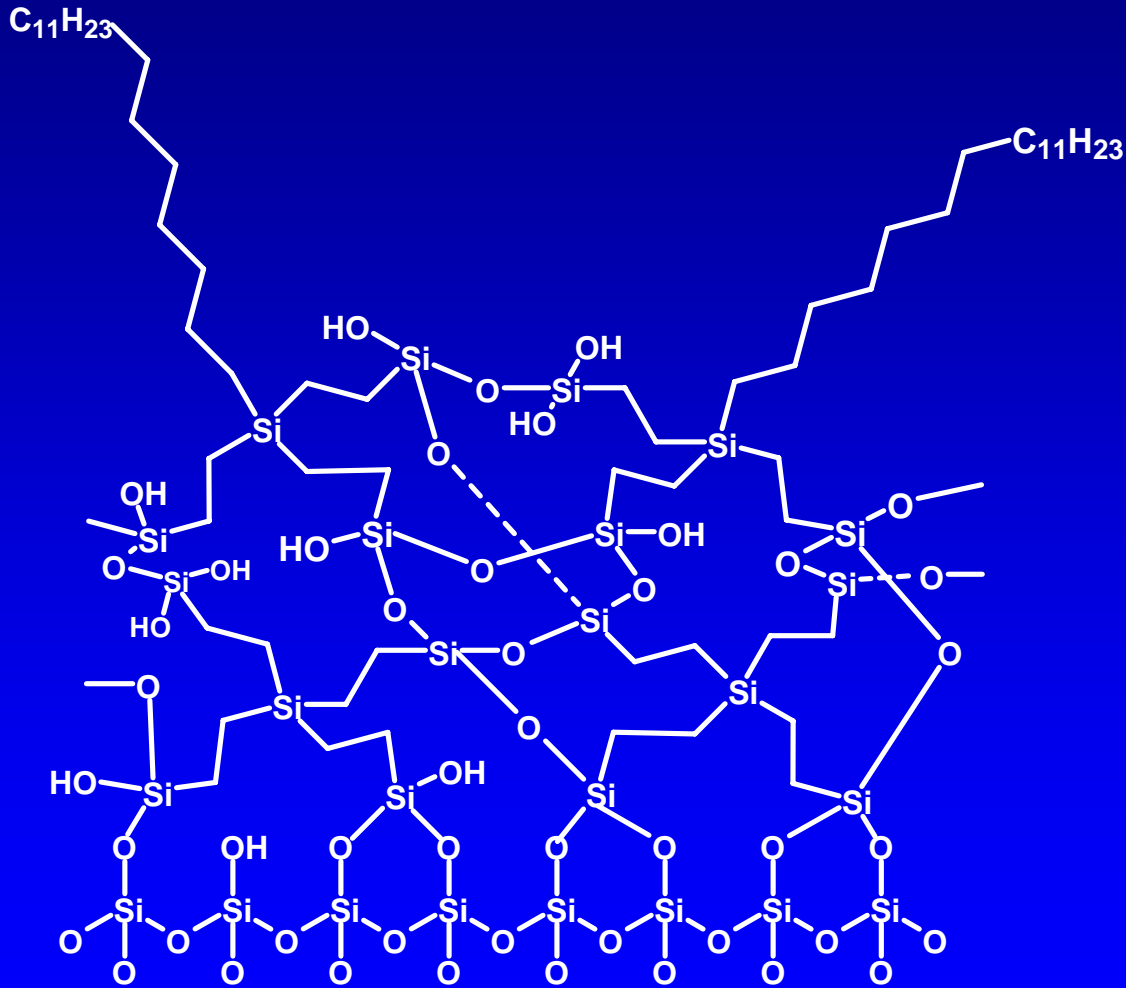

US and International Patents Pending

SFC Test of pHidely™C₁₈ Column

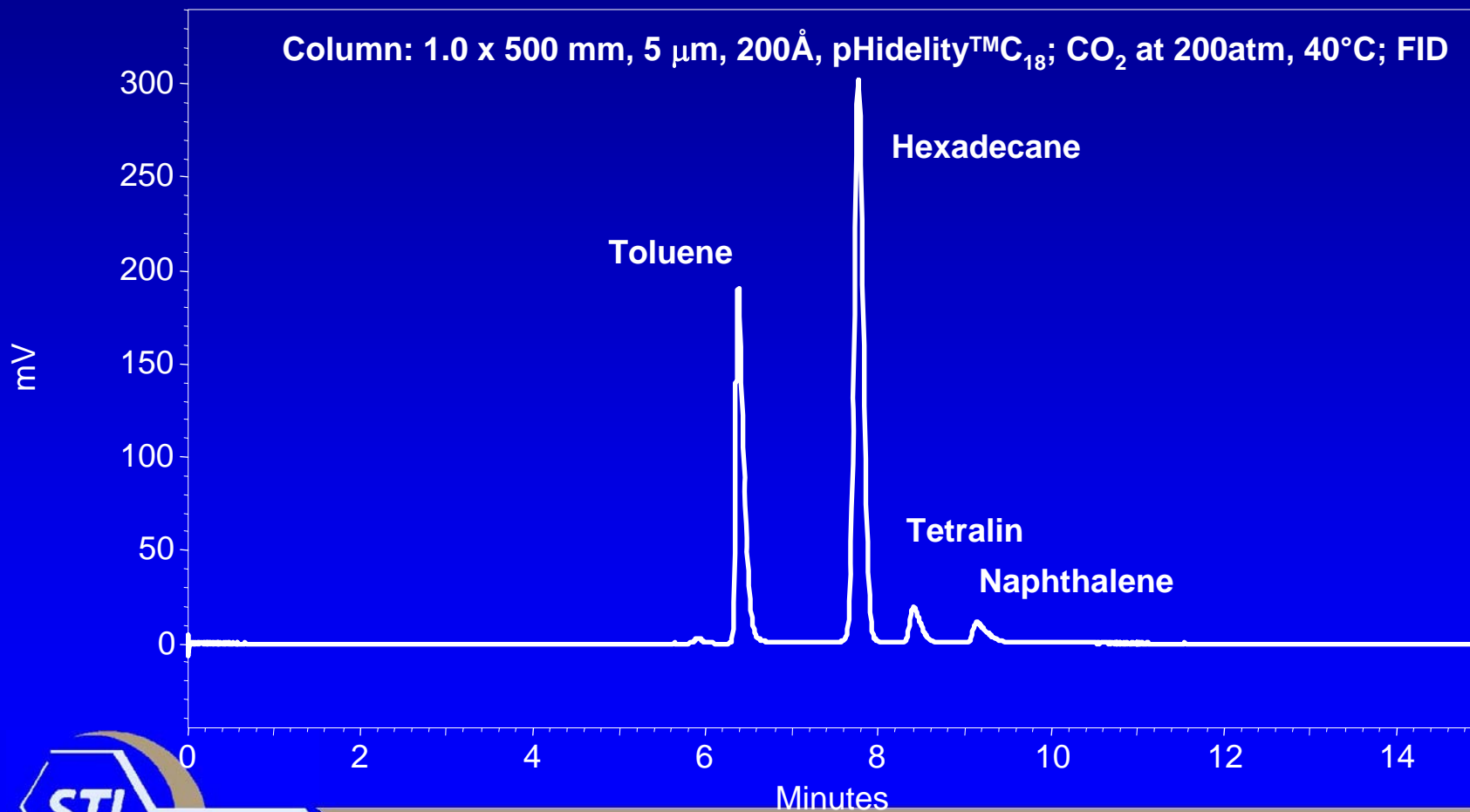


pHidely is a Trademark of Restek Corporation

Selerity Polaratherm™ Series 9000 Total Temperature Controller

- Used in the LC portion of this study
- Forced air oven and chiller
- Isothermal and thermal gradient operation
 - Sub-zero to 200°C
 - Thermal gradients up to 30°C/min
- Mobile phase preheating and pre-cooling
- Peltier effluent temperature control
- Vapor sensor
- Compatible with any HPLC system



Resolution Test

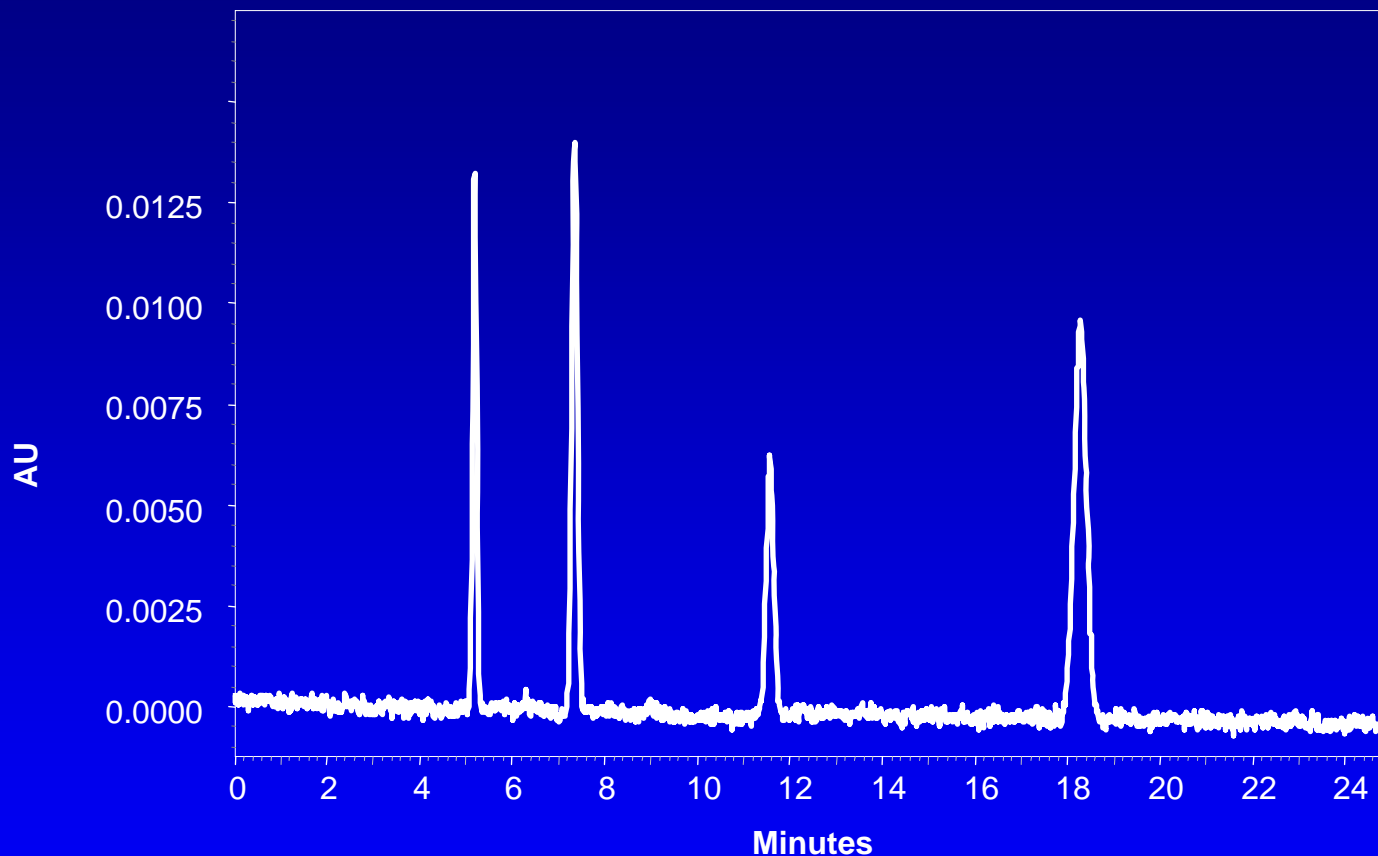


Column: 1.0 x 500 mm, 5 μ m, 200Å, pHidelity™ particles, 50°C
UV 254nm, 170nL cell; 60nL injection; Flow: 0.05 ml/min 50%ACN

Elution: uracil, phenol, benzene, naphthalene; Reduced plate height: ~4




Pressure Drop vs. Temperature

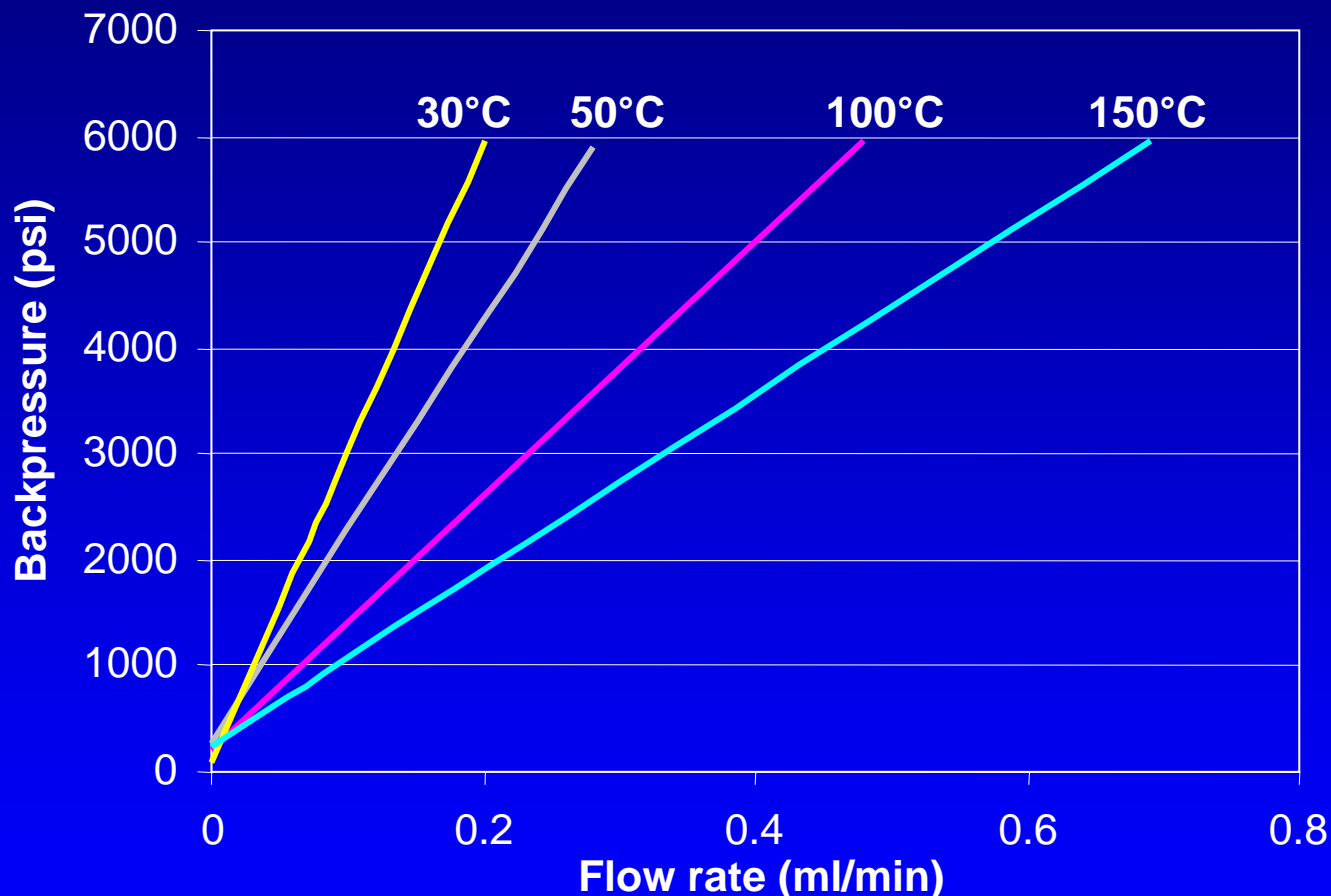

50%ACN 1x500mm pHidely™/Blaze™ column 5um particles



High Speed Separation with a 1 x 500 mm Column

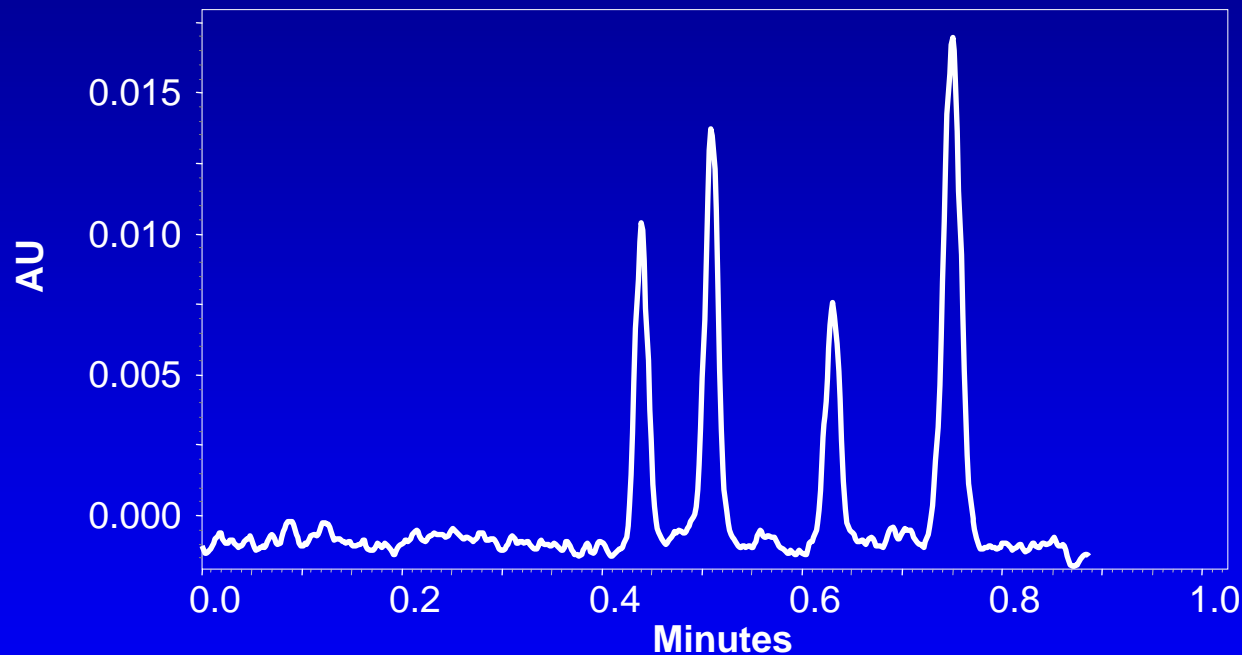


Column: 1.0 x 500 mm, 5 μm , 200 \AA , pHidelity™ particles, 150°C

Detection: UV 254 nm, 170nL cell, Flow: 0.4 ml/min 50%ACN

Elution: uracil, phenol, benzene, naphthalene



Sleep Aid Separation

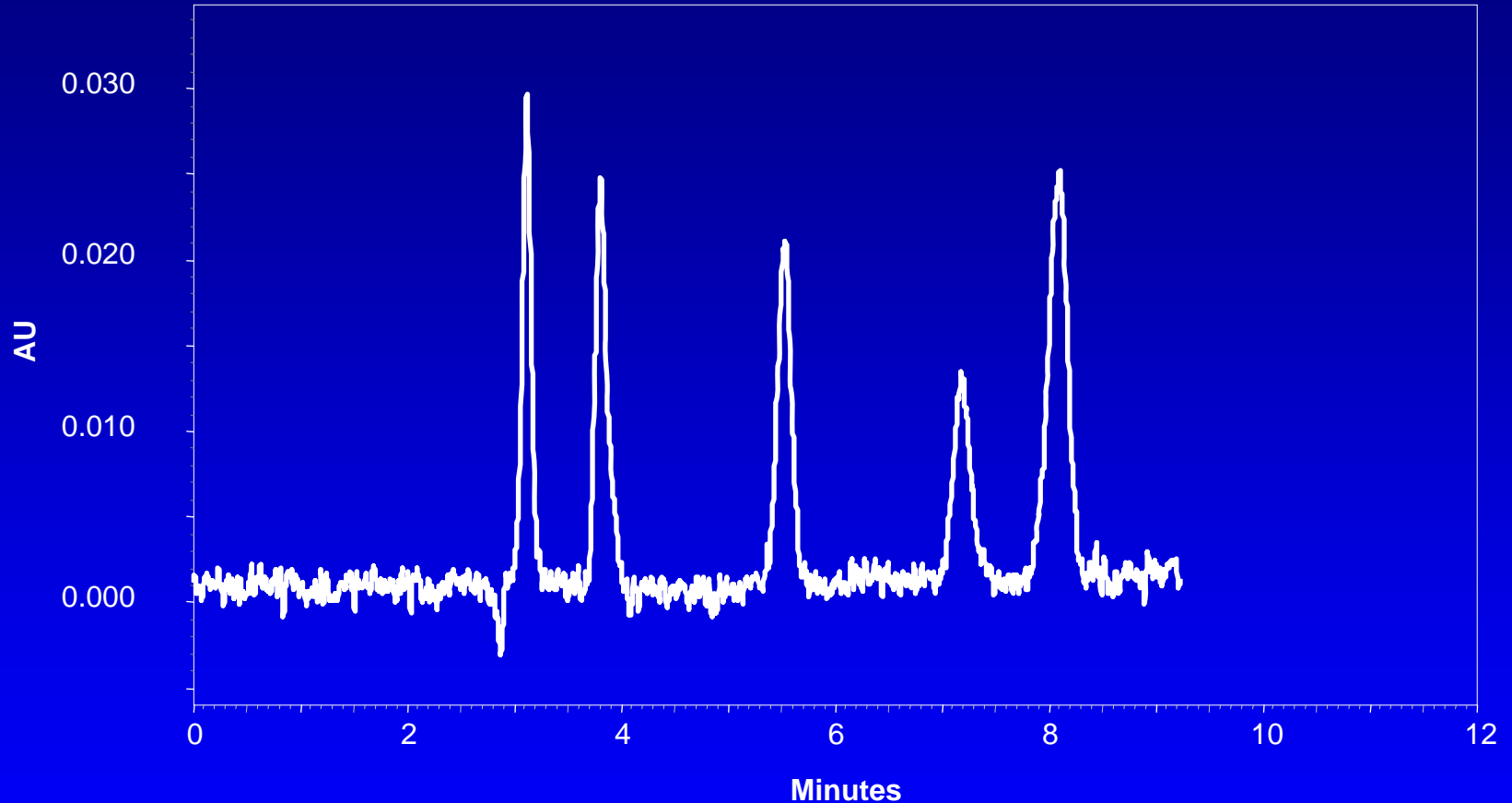


Column: 1.0 x 500 mm, 5 μ m, 200 \AA , pHidalityTMC₁₈ particles, 50 $^{\circ}$ C; UV 254 nm
170nL cell, 60nL Injection; Flow: 0.1 ml/min 35%ACN, 0.2%TFA

Elution: doxylamine, LunestaTM, AmbienTM, diphenhydramine, SonataTM



Analgesic Separation

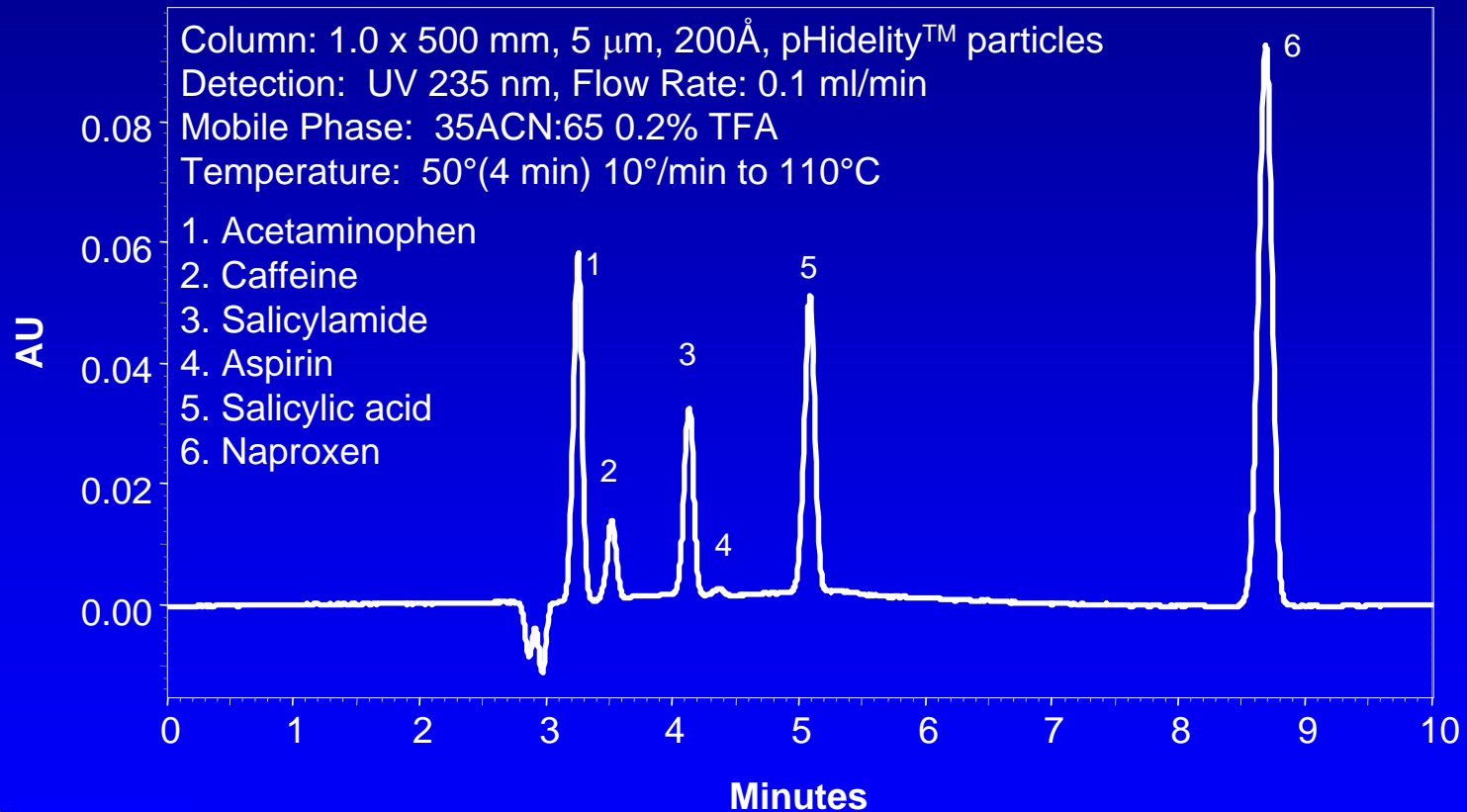

Conclusions

- Polished stainless steel tubing is capable of producing narrow-bore columns in long lengths with good efficiency.
- Stable narrow-bore columns for reversed-phase chromatography can be used with wider pH and temperature ranges.



 **Turn up the Heat!** 

www.selerity.com

Partnering with Restek Corporation

