

Optimized Instrument Parameters for a High Temperature HPLC System with Flame Ionization Detection

Jody Clark, Stephanie J. Marin, Brian A. Jones and
W. Dale Felix

Selerity Technologies, Inc.
2484 W. Custer Road
Salt Lake City, UT 84104



www.selerity.com


High Temperature Liquid Chromatography (HTLC) Advantages

The major advantages are

- Decrease analysis time (increase linear velocity)
- Increase efficiencies and resolution
- Tune selectivity with temperature
- Decrease organic solvent usage



Elution Strength Illustration: Separation of Steroids Using Water as the Mobile Phase

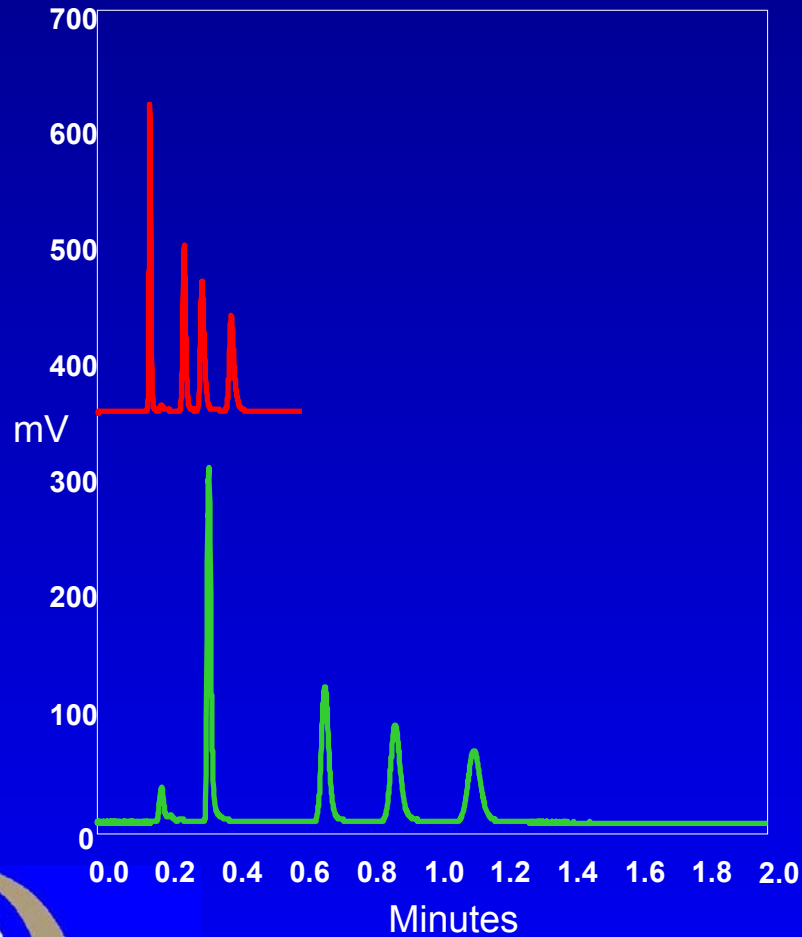

Column: ZirChrom PBD, 3 μ m
100 X 4.6 mm

Detection: UV 254 nm

Flow Rate: 6.0 mL/min

Mobile Phase: Water

Temperature: 200°C

Flow Rate: 3.0 mL/min

Mobile Phase: 25:75 acetonitrile:water

Temperature: 50°C

Elution Order:

Uracil


Androstadienedione

Androstenedione

Epitestosterone



Solvent Polarity as a Function of Temperature

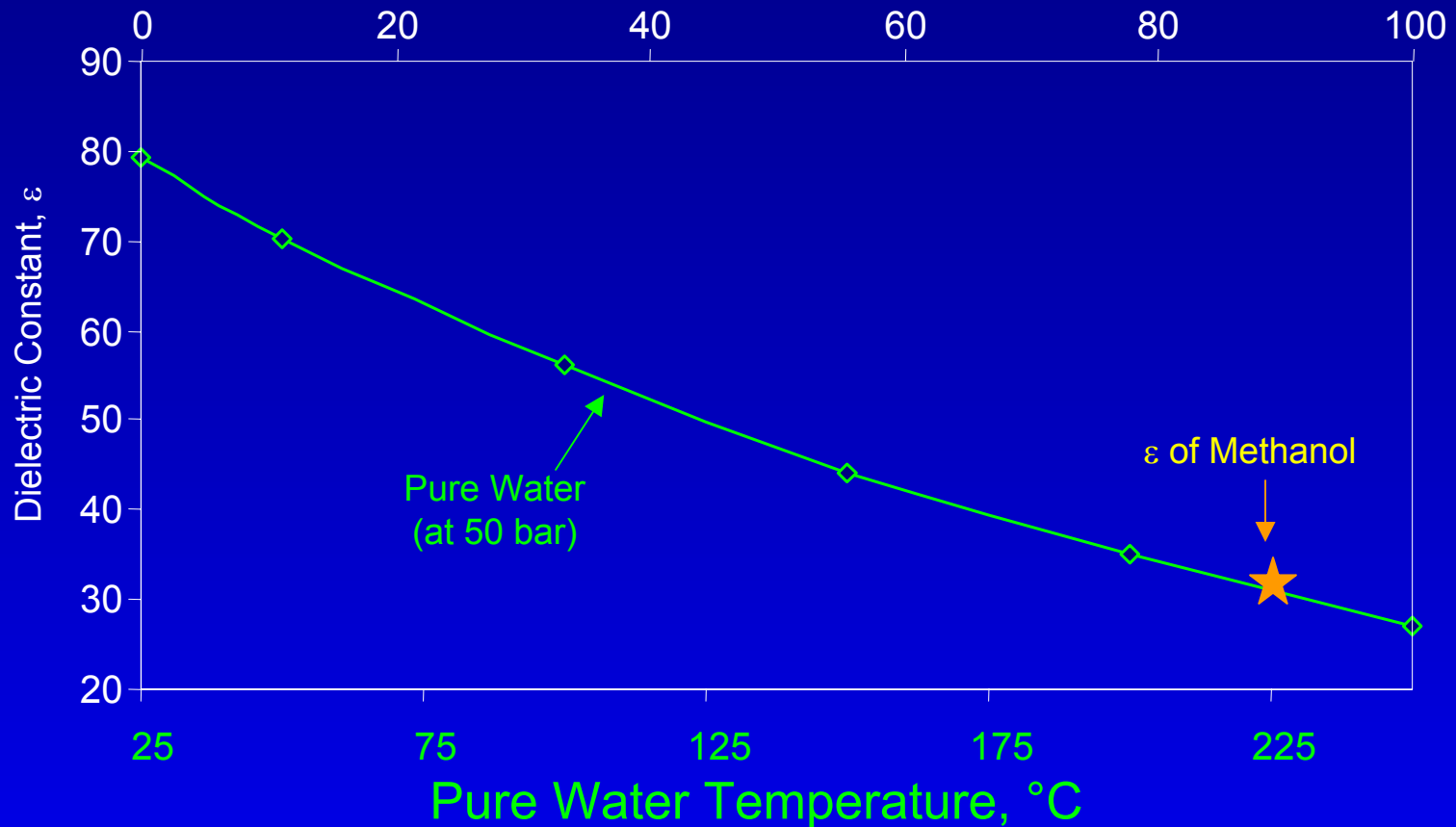

Data from Y. Yang et al. *J. Chromatogr. A* **810** (1998) 149.




Why Use Water as a Mobile Phase?

- Inexpensive
- Readily available
- Non-polluting
- Transparent to most detectors including UV, FID, and NMR (D_2O)
- Elution strength can be controlled by varying temperature



Temperature Affects of Hydrogen Bonding

- Increasing temperature-
 - Increases intermolecular distance
 - Weakens hydrogen bonds

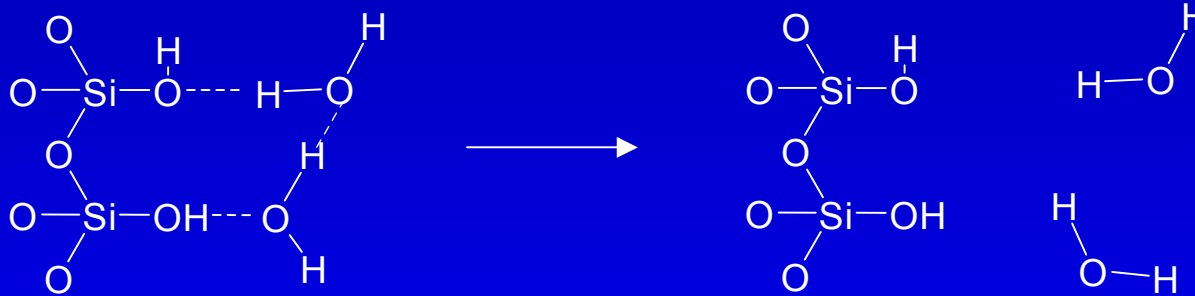

Advantages of Using Flame Ionization Detection?

- Mass flow sensitive detector
- Response is proportional to the number of carbon atoms being burned
- Wide linear range, about 10^8
- Responds with high sensitivity to organic compounds




Instrumentation

- Aquachrom temperature programmable HPLC oven with FID
- Mobile phase preheater
- Split assembly
- Mobile phase pure water
- Eldex syringe pump
- Hypercarb™ column with stainless steel end fittings



FID Design

- Standard GC type design with grounded jet
- 0.020" internal diameter flame jet
- Temperature limit 400°C

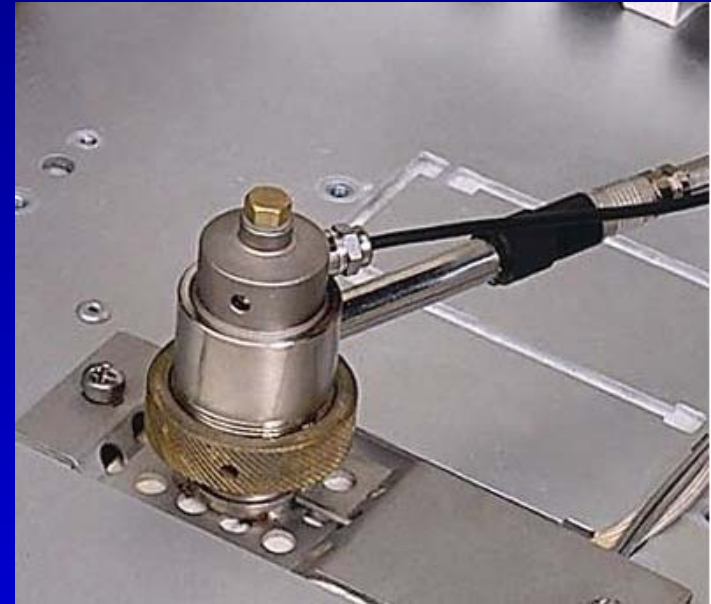


Optimization Parameters

- Detector interface
- FID conditions
- LC column sizes
- Pump flow rate




Restrictor Types for Superheated Water

- Stainless steel (silica degrades)
- Create backpressure to keep water in liquid form
- Pressure drop short distance (heated zone)
- Introduce water into the flame in aerosol

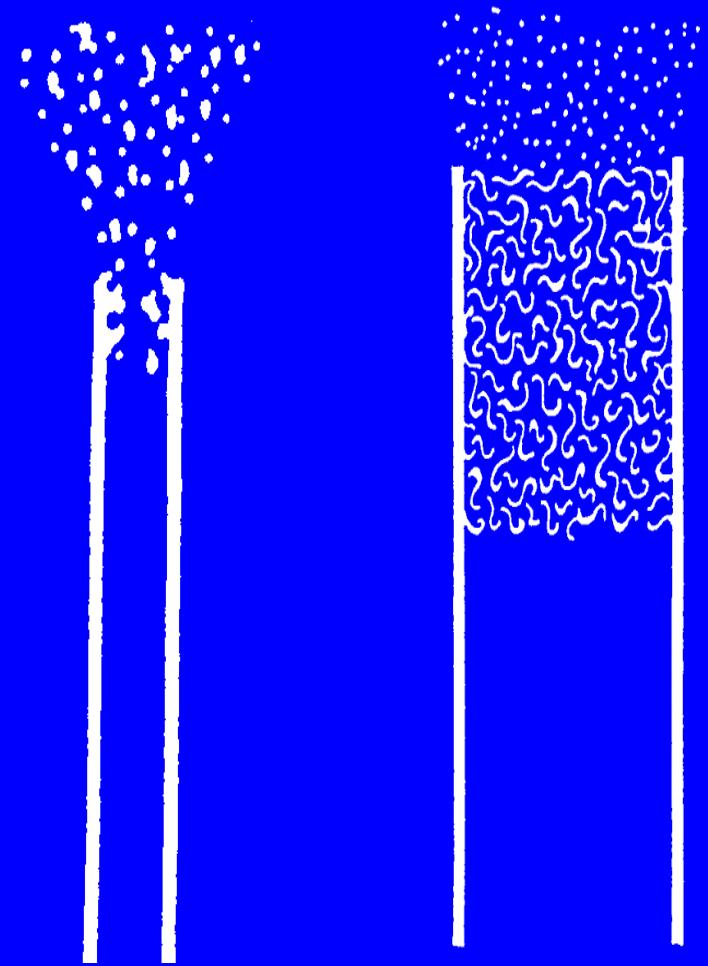



Selerity Restrictor

- Introduce water into flame in aerosol form

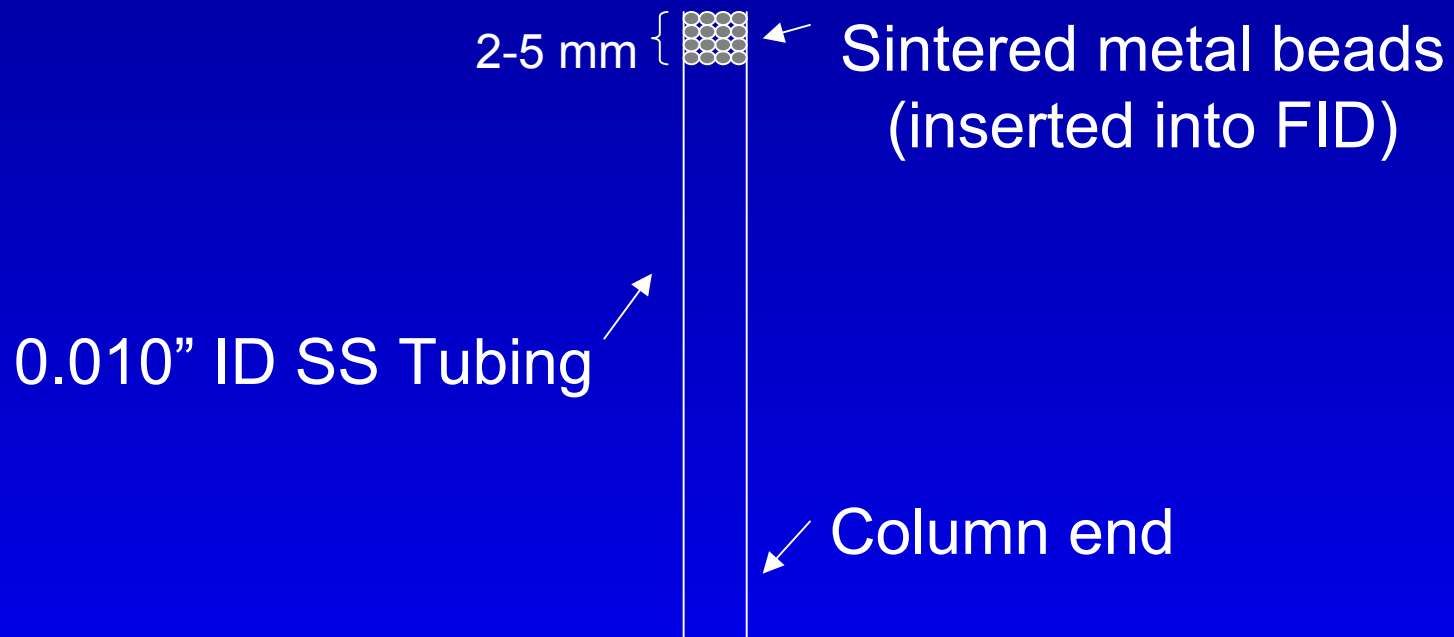

Linear vs. Metal Restrictor

- Direct into flame jet (no column)
- Methanol
- Temperature ramp: 50-200°C @ 20°C/min.

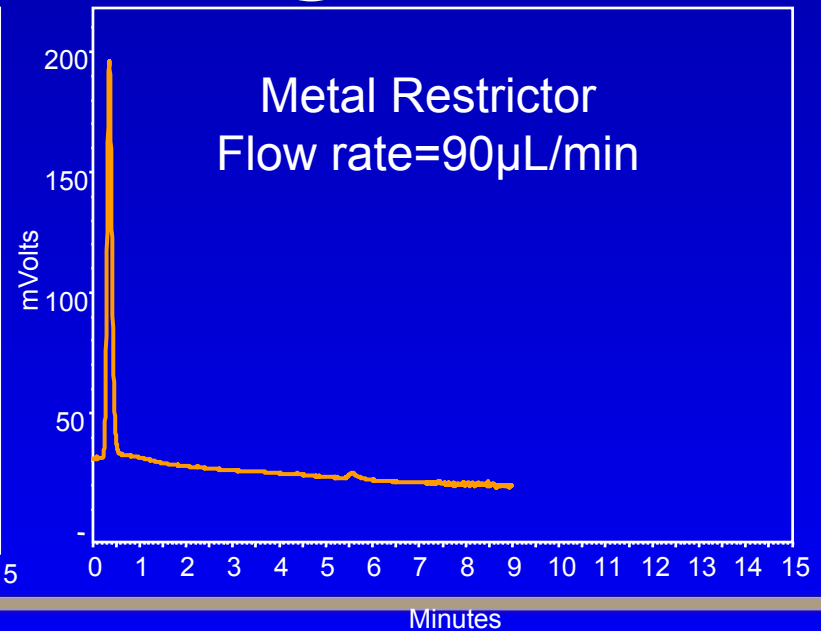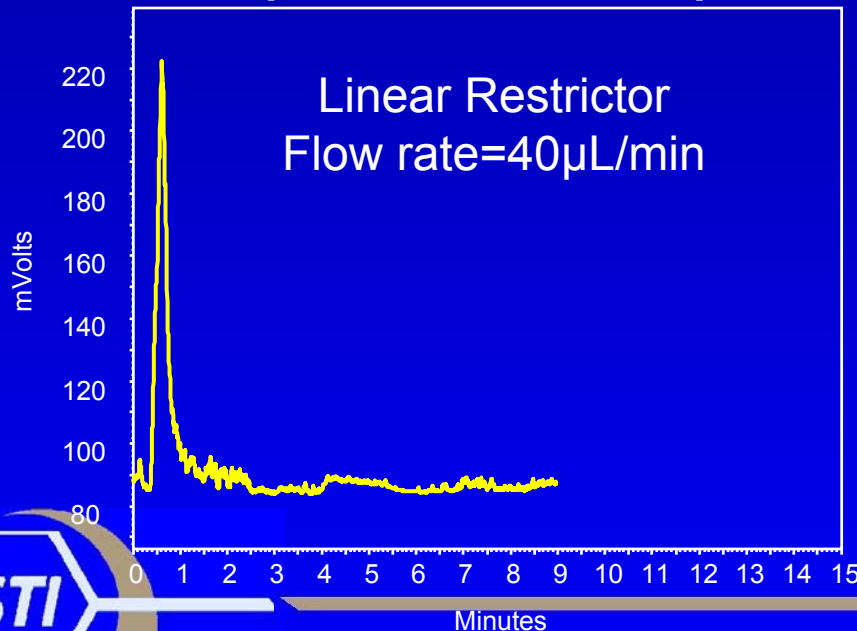

Factors Affecting Sensitivity

- Volume of water and sample introduced into FID
- FID temperature
- Restrictor position
- Detector gases



Superheated Water Flow into Flame

- Started with SFC-FID conditions
- Direct connection from injector to FID (eliminate column interference)
- Oven program from 50 to 200°C

Results

- Maximum amount of water flow equals 100 $\mu\text{L}/\text{min}$




FID Temperature


- FID temperature must be above 375°C
- No significant change in sensitivity or peak shape at FID temperatures of 375 and 400°C



Baseline Difference Between 0.5 cm and 1.0 cm

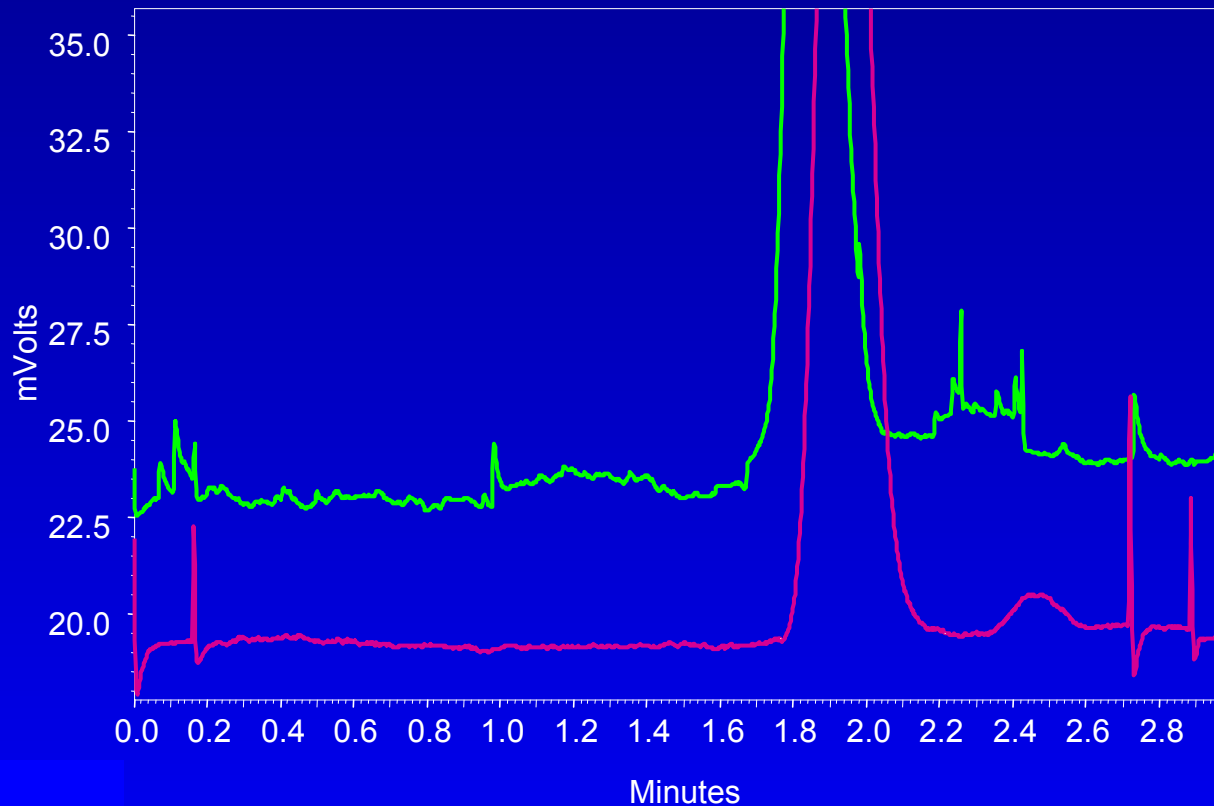

Distance of Restrictor From Tip of Flame Jet

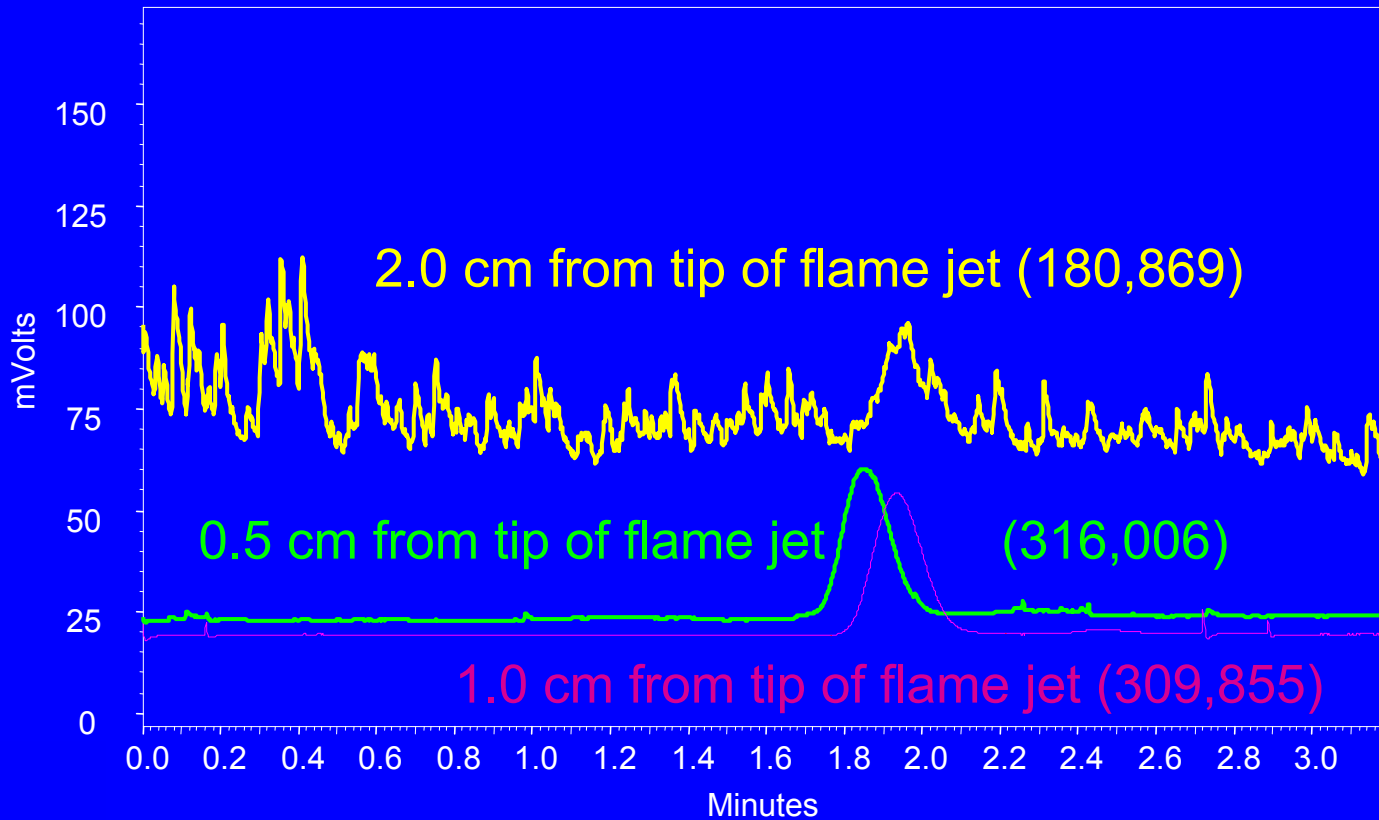


Detector Gas Flow Rate (H₂)


- Oven program 50 to 200°C

H ₂ flow rate	Area	Variance
75 mL/min	316,006	Reference
100 mL/min	283,637	10 %
115 mL/min	237,095	25 %

100 mL/min

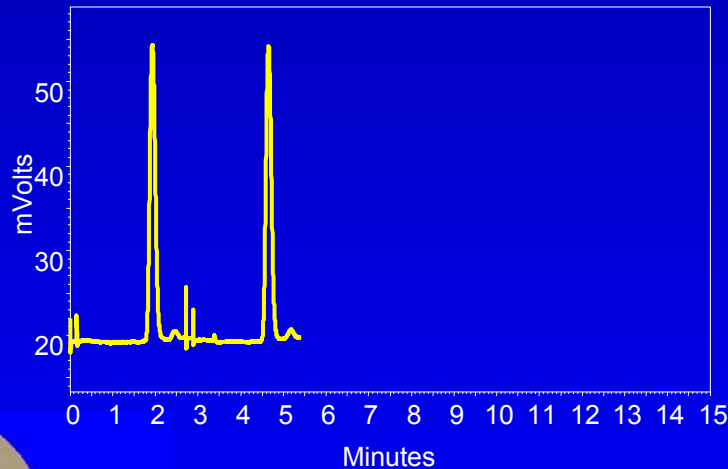


115 mL/min

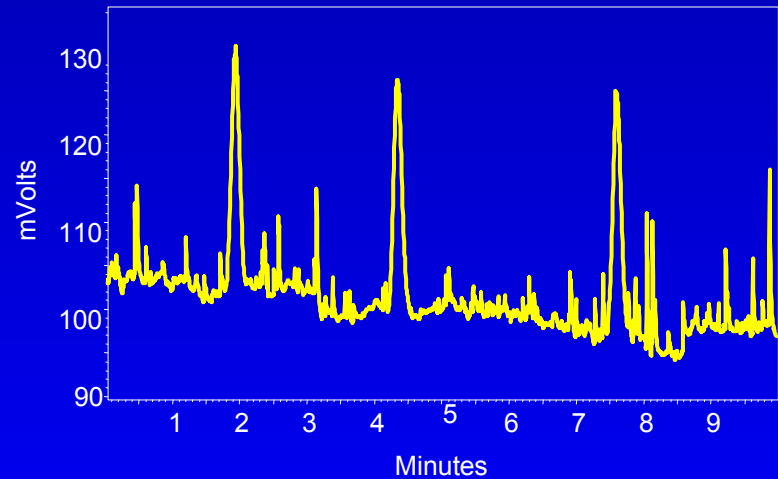


Effect of Make-up Gas on Sensitivity (Nitrogen)

- 8.4% loss of area without nitrogen

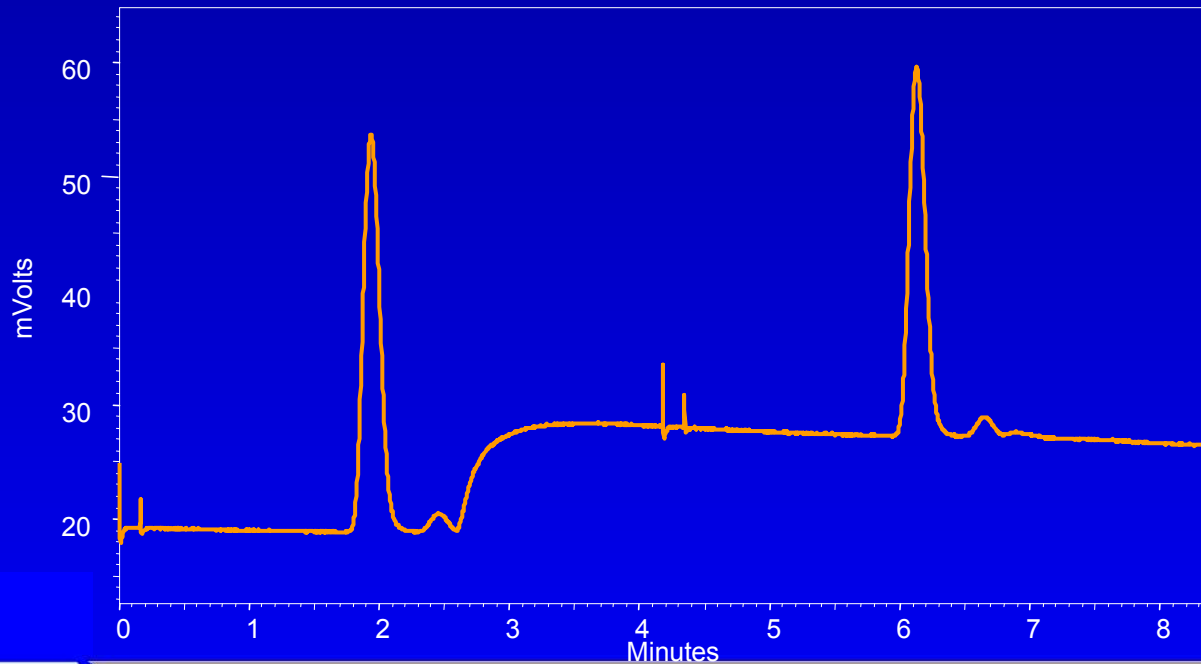


Sensitivity

- 10.4 ng of methanol on column (Area=28,302)
- Optimum conditions

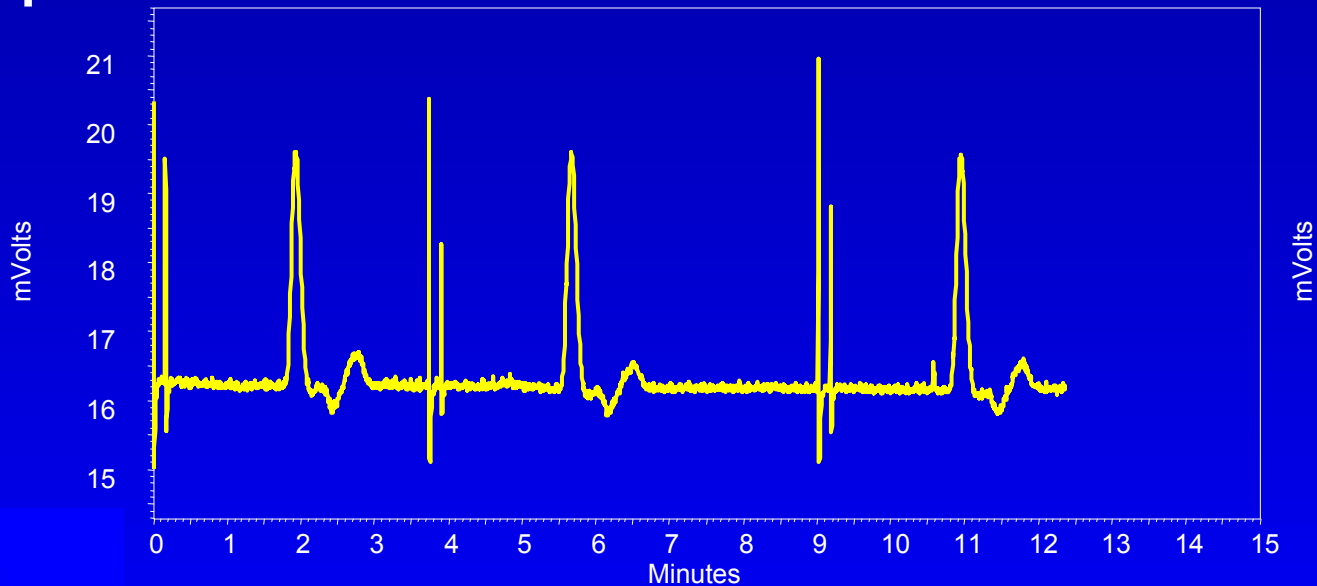

Detector Linearity

Concentration Range= 0.00208-0.415 mg/mL

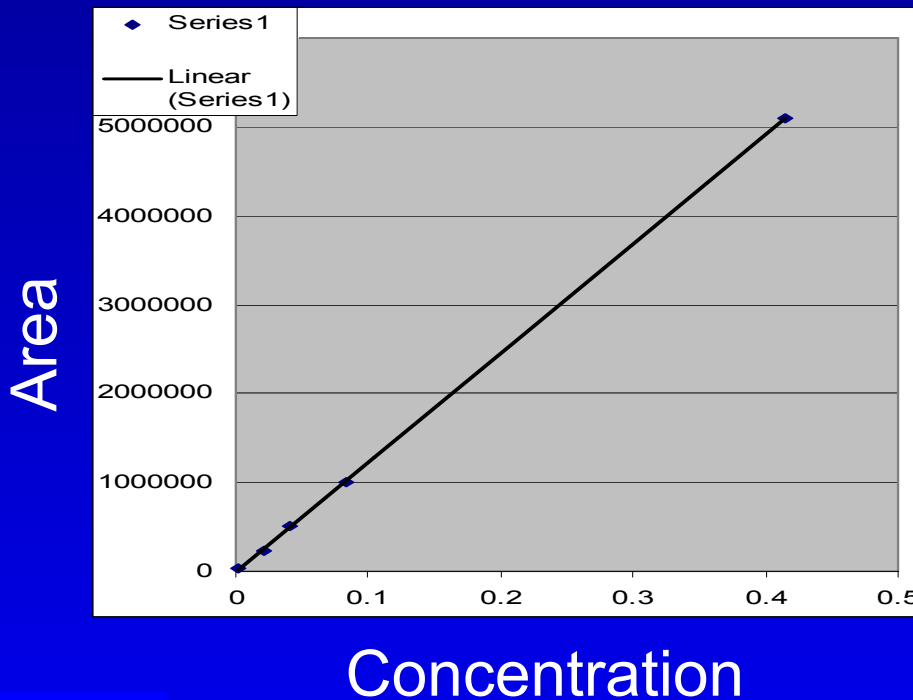

Regression Coefficient

$$R^2=0.99996159$$



Traditional LC Column Sizes and Optimal Pump Flow Rates

- 1.0 mm ID
 - Optimum flow = 50 $\mu\text{L}/\text{min}$
 - Superheated water into flame 50-100 $\mu\text{L}/\text{min}$
- 2.1 mm ID
 - Optimum flow = 200 $\mu\text{L}/\text{min}$ (must split flow)
 - Up to 1.0 mL/min with split
- 4.6 mm ID
 - Optimum flow = 1.0 mL/min
 - 10% sample elutes to FID
 - Use concentrated samples



Split Assembly

- 40 μm ID fused silica




Flow Rate vs. Area (2.1mm ID Columns)

Water flow rate	Area	Theoretical split	Length split tubing(40 µm ID)
200 µL/min	2,880,393	50%	25 cm
300 µL/min	1,523,720	67%	20 cm
400 µL/min	1,381,774	75%	10 cm
500 µL/min	1,274,604	80%	10 cm
600 µL/min	1,056,077	83%	10 cm
700 µL/min	938,846	86%	5 cm
800 µL/min	648,870	88%	5 cm
900 µL/min	560,189	89%	5 cm
1000 µL/min	562,955	90%	5 cm




1 mg/mL Methanol

Glycols on 1.0mm ID Hypercarb™

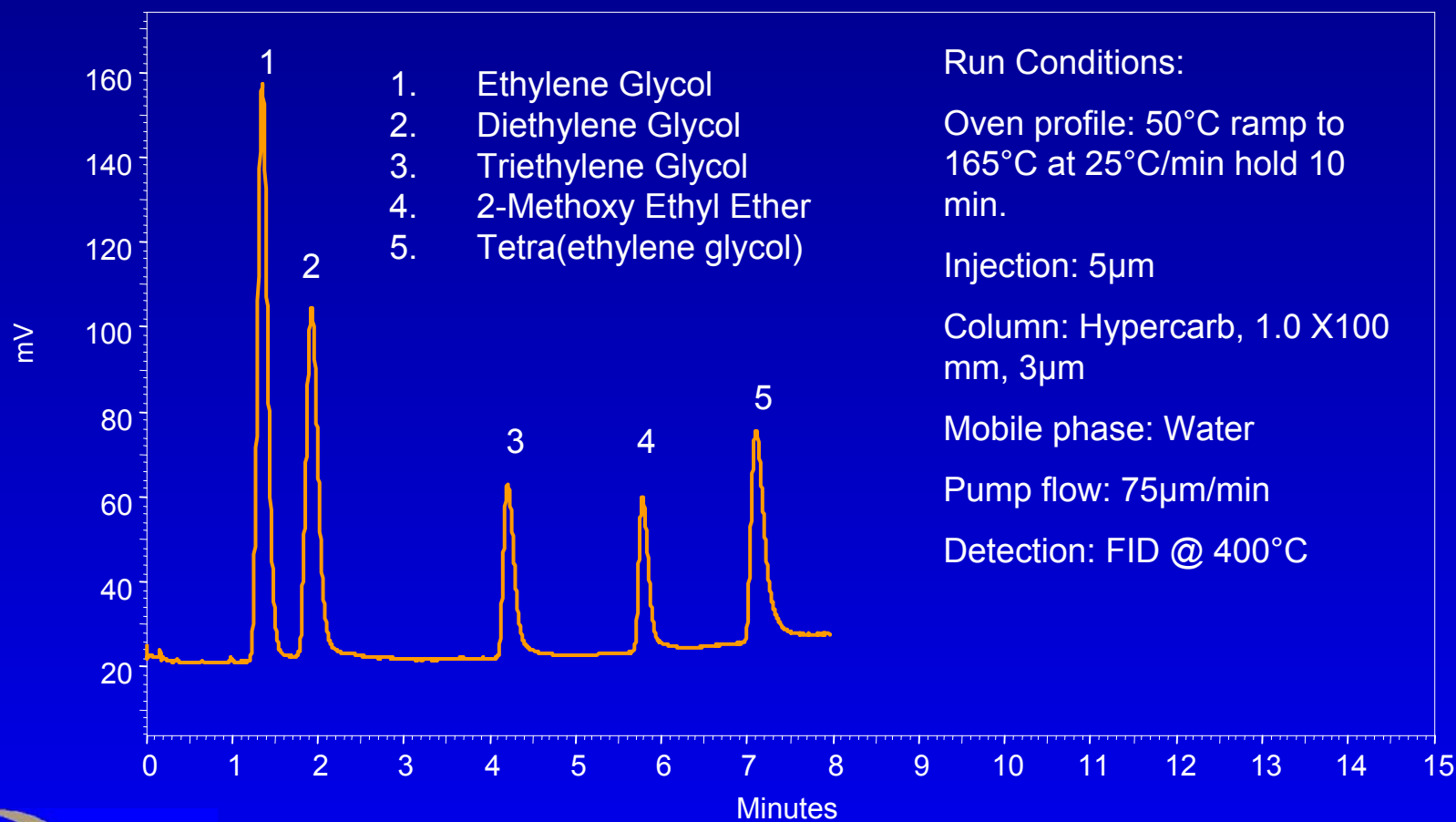

Alcohols on a 2.1mm ID Polymer Column

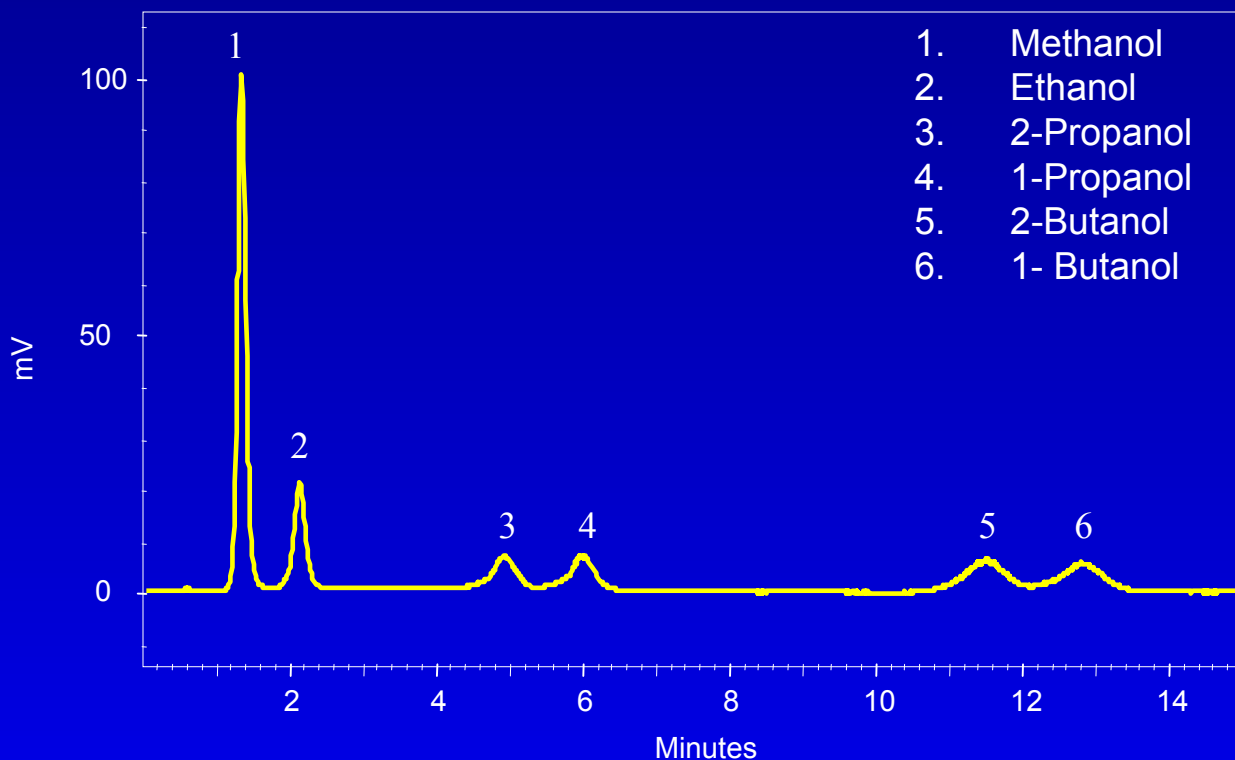

Run Conditions:

Oven profile: 50°C ramp to 135°C at 10°C/min hold 10 min.

Column: Jordi DVB, 2.1 X100 mm, 5µm

Mobile phase: 0.025%TFA

Split flow: 40µm ID fused silica back pressure

Pump flow: 0.500 mL/min

Detection: FID @ 400°C



Conclusion

- Optimum conditions depend on separation goals
- Sensitivity: use low hydrogen flow, low pump flow rates and 1.0 mm columns
- Fast analysis or dual detection: use higher hydrogen flow rates, higher pump flow rates, larger columns and split flow



Acknowledgements

- Thermo Electron Corporation
- Nelu Grinberg, Ph.D., Boehringer Ingelheim Pharmaceuticals, Inc.

